

Achievement of 1 Mcps Output Rate from a Large Area Silicon Drift Detector

L. Feng, V.D. Saveliev, M. Takahashi, S. Barkan, E.V. Damron, C.R. Tull

SII NanoTechnology USA, Northridge, CA 91324

We are reporting here for the first time the achievement of one million counts per second (1 Mcps) output count rate (OCR) from our large area (~50 mm²) single element silicon drift detectors (the Vortex[®]) using a commercially available digital pulse processor, the Mercury Digital X-ray Processor from XIA (X-ray Instrumentation Associates, LLC).

The advantages and high performance of the Vortex[®] silicon drift detectors have been extensively reported. However, their highest achievable OCR has been limited to 500 - 600 Kcps in the past, due to limitations of the previous generation of pulse processors. Old processors offer a minimum processing peaking time of only 0.25 μ s, with a limited capability in pile-up rejection (PUR). This made it impractical to break through the 500-600 Kcps OCR limit, although the Vortex[®] detectors feature a very short rise time (< 100 ns) which is suitable for a much shorter processing peaking time. Recent advances in digital pulse processing technology has led to a new generation of pulse processors which typically run at much higher clock speed (\geq 50 MHz) and offer much shorter processing peaking times (\leq 0.1 μ s) with improved PUR performance, making it possible to exceed the previous OCR limit. Using XIA's Mercury processor, for example, we achieved 1 Mcps output rate with 0.1 μ s processing peaking time. This is a significant improvement, especially for applications where OCR performance is critical, such as in fast X-ray mapping, ultra-trace elemental analysis, EXAFS, etc. Figure 1 shows a series of Mn K spectra acquired for 0.1 μ s peaking time at various OCR from a Mn sample excited with a Rh X-ray tube using 40 kV voltage. For up to ~3 Mcps input count rate (ICR), both peak position and resolution are reasonably stable: the peak position change is less than 1% while the Mn K _{α} resolution varies within the range of 345 - 386 eV. The acquisitions were set to have a moderate PUR performance as exhibited in Figure 1.

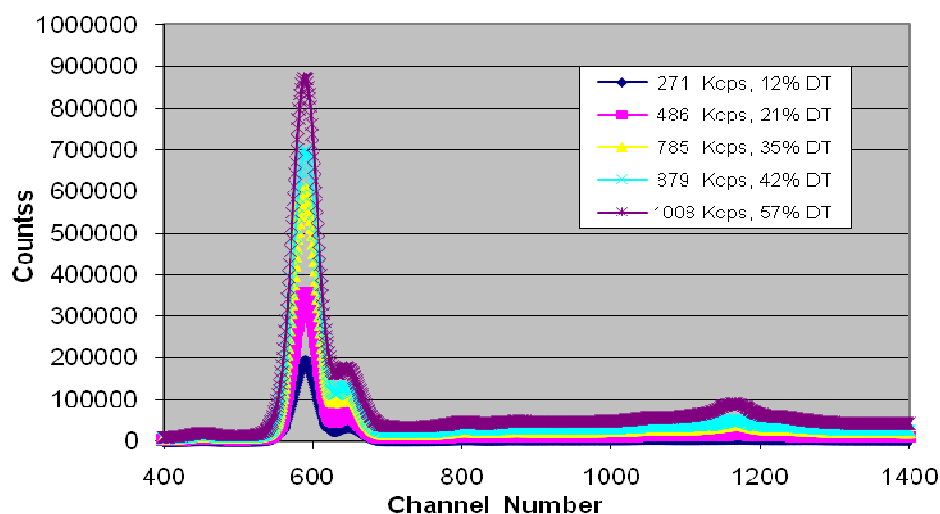


Figure 1. Mn spectra acquired using 0.1 μ s peaking time at various OCR